



Oxygraph-2k Manual


Mitochondrial Physiology Network 19.18(D01): 39-48 (2014)
D: www.bioblast.at/index.php/MiPNet19.18D_O2k-Calibration

©2014 OROBOROS®
Version D01: 2014-10-01


O2k calibration by DatLab

Gnaiger E


OROBOROS INSTRUMENTS Corp
high-resolution respirometry
Schöpfstr 18, A-6020 Innsbruck, Austria
Email: erich.gnaiger@oroboros.at
www.oroboros.at

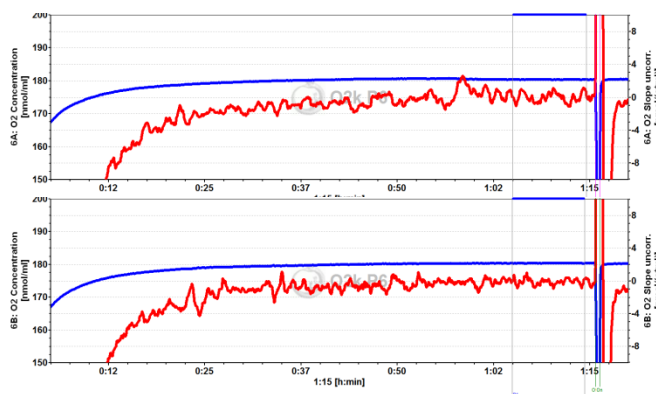

OROBOROS Oxygraph-2k

Section		Page
	1. Experimental oxygen calibration	40
	2. DatLab-calibration of oxygen sensors	42
	3. MultiSensor calibration	46

Summary Accurate calibration of the oxygen sensor depends on (1) incubation medium equilibrated at oxygen partial pressure of air, achieved in the Oxygraph-2k (O2k) at defined temperature, continuously recorded total gas pressure (barometric pressure), and thermodynamic equilibrium between the gas and aqueous phase; (2) zero oxygen calibration; (3) high stability of the signal of the polarographic oxygen sensor (POS), tested for sufficiently long periods of time; (4) linearity of signal output with oxygen pressure, achieved with the POS in the range between oxygen saturation and zero oxygen pressure; and (5) accurate oxygen solubilities for aqueous solutions for the conversion of partial oxygen pressure into oxygen concentration ([MiPNet06.03](#) with O2k-SOPs). The standard oxygen calibration procedure is described for O2k high-resolution respirometry with the automatic calibration routine by DatLab. MultiSensor calibration by DatLab is explained briefly.

 Before completing the Oxygen calibration [F5] in DatLab, oxygen concentration [$\mu\text{M}=\text{nmol/ml}$] is displayed on the basis of default calibration settings. Oxygen flux [$\text{pmol}\cdot\text{s}^{-1}\cdot\text{ml}^{-1}$] is accurate in terms of high-resolution only after oxygen calibration and instrumental background calibration ([MiPNet19.18E](#)). Default

calibration values are replaced by experimental calibration, whenever sufficient stability of the calibration cannot be assumed, or when previous calibration conditions do not apply. The recorded raw signal, R_t , is converted to oxygen concentration, $c_{O_2,t}$ [μM], or partial pressure, $p_{O_2,t}$ [kPa or mmHg]. In DatLab, calibration parameters can be edited any time during the experiment.

1. Experimental oxygen calibration

The polarographic oxygen sensors (OroboPOS, POS) are calibrated by a two-point calibration, routinely achieved at air saturation and zero oxygen concentration. Accordingly, static calibration involves the determination of the constant signal of the POS recorded at 100% and 0% air saturation (R_1 and R_0) under the particular experimental conditions (temperature, signal amplification by electronic gain, polarization voltage, stirring speed, medium).

1.1. Air calibration

O2k-SOP

Air saturation is achieved by stirring the aqueous medium in contact with air in the O2k-Chamber without sample, following the procedures below and in [MiPNet06.03](#).

1. Add incubation medium into the chambers, with an excess volume of at least 0.1 ml above the experimental chamber volume (2 ml), to fill the O2k-chamber and injection capillary of the stopper when it is fully inserted (closed). The volume does not have to be accurate, as long as it is above the minimum volume. Switch on the stirrers either during or after addition of the medium.



Stopper-spacers used for air calibration with white PVDF stopper (Chamber A) or black PEEK stopper (Chamber B).


2. Insert the stoppers slowly to their volume-calibrated position ([MiPNet19.18A](#)). Siphon off excess medium ejected through the injection capillary and collected in the well of the stopper. Then lift the stoppers using the stopper-spacer tool, leaving a gas volume above the liquid phase for final air equilibration.

The central level of the gas phase remains above the rotating stirrer bar, preventing bubbles and foam from being formed which would block gas exchange. To

ensure a well defined p_{O_2} in the gas phase, the gas volume has to be renewed (exchanged for air), if the medium was originally not near air saturation. This is achieved simply by fully inserting and re-opening the stopper. Equilibration is a slow process: stability should be reached within one hour (figure below). A stirrer test [F9] can be performed during equilibration ([MiPNet06.03](#)).

3. After stabilization of the POS signal, the recorded signal at air saturation, R_1 , is about 2-3 V at Gain 1. A signal of 1 V corresponds to a signal current of the POS of 1 μA (corresponding to 2 V at Gain 2). Under all experimental conditions, the raw signal must be <10 V.

Continue recording for 3-10 min to check for signal stability. You may proceed at this point with an O2k-background test ([MiPNet14.06](#)).

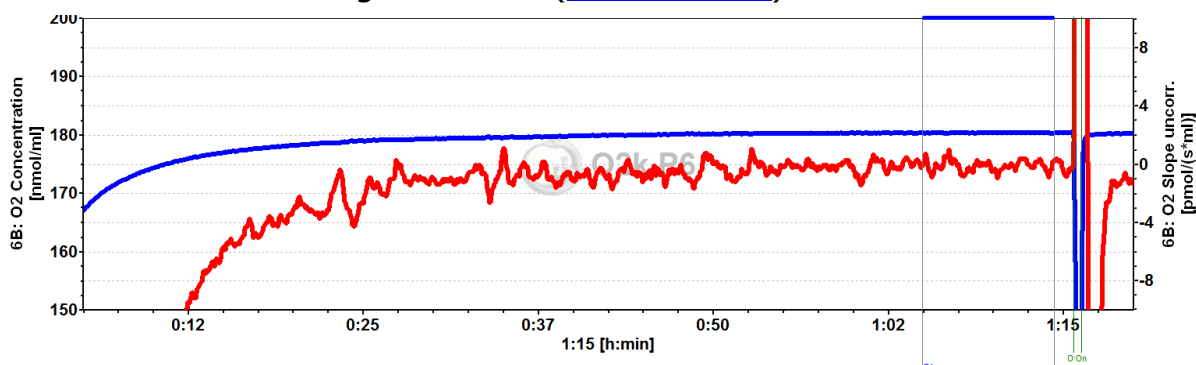

Oxygen concentration (blue plot; 180 μM ; full scale 50 nmol/ml or 50 μM) over 1 h after switching on the Oxygraph-2k (Power-O2k P6, chamber B) and setting the experimental temperature at 37 $^{\circ}\text{C}$, with medium stirred for equilibration with a gas phase of air at 575 m altitude. The **red plot** is the negative slope of oxygen concentration over time [$\text{pmol}\cdot\text{s}^{-1}\cdot\text{ml}^{-1}$] on the right Y-axis, with zero in the middle position. A slope of zero (for 'O2 Slope uncorrected') indicates a constant O_2 signal over time. 2014-02-19 P6-01.DLD

1.2. Zero oxygen calibration

4. Zero oxygen calibration is achieved best with mitochondria or cell suspensions by allowing complete oxygen depletion. Alternatively, titrate zero solution (Na-dithionite, [OroboPOS-Service Kit](#)) into the chamber, which is part of the automatic TIP2k-supported instrumental background test ([MiPNet14.06](#)). The stable zero signal, R_0 , should be $<2\%$ of the signal at air saturation, but $<5\%$ is acceptable. Most importantly the zero signal must be stable.



go Bioblast » [MiPNet06.03 POS-Calibration-SOP](#)

2. DatLab-calibration of oxygen sensors

2.1. Graph layout for calibration

Graph Layout: **"O1 Calibration Exp. Gr3-Temp"**

Calibration experiment with temperature and Peltier power in Graph 3.

This is typically the first layout used after switching on the O2k. Oxygen concentration (blue lines, left Y-axis) and O2 slope uncorrected (red lines, right Y-axis) are displayed on the top graph for the left chamber, and below for the right chamber. The third graph (bottom) shows the block temperature on the left Y axis and the Peltier power on the right Y axis. Only when both temperature and Peltier power are constant, the chambers have reached thermal equilibrium. The next step is to observe equilibration of the oxygen signal with a defined gas phase above the stirred aqueous phase ('open' chamber; usually with air as the first step) to perform an oxygen calibration.

If anything unusual is observed (always zero flux, jumping signals), the layout "Z Trouble Shooting" should be used.

Graph Layout: **"O2 Background Experiment"**

For recording O2 sensor calibration and a test for instrumental background oxygen flux.


For each chamber, 'O2 Concentration' and 'O2 Slope uncorrected' are displayed on the left and right Y-axis, respectively. 'O2 Slope uncorrected' is the negative slope of oxygen concentration (multiplied by 1000 to convert to units [pmol/ml]) over time [s]. No correction is applied for instrumental background oxygen flux. Zero flux in the 'open' chamber at air calibration indicates stability of the oxygen signal. After closing the chamber, 'O2 Slope uncorrected' deviates from zero as a function of the oxygen consumption of the polarographic oxygen sensor and of oxygen diffusion into or out of the chamber.

2.2. Mark

Mark a section of the experiment at air saturation, when signal stability is reached. This should be done real-time to save default calibration information. Corrections are possible after disconnecting from the O2k. For calibration, follow steps (1) to (7) illustrated on the following graph:


1. Select a graph by a L[⌘] click (left mouse button) into the graph, or directly by step 2.
2. Select the oxygen signal as the active plot by a L[⌘] click on Y₁ in the figure legend on the right of the graph. The active plot is highlighted.
3. Only if 'Mouse Control: Zoom' mode has been activated: Select "Mouse Control: Mark" in the Graph menu, or press [Ctrl+M].
4. Set a mark [Shift+L[⌘]]: Hold [Shift] and click the left mouse button, move the cursor along the time axis, release the left mouse button at the end of the section to be marked.

Remove a section of the mark or the total mark [Shift+R[⌘]]: Move the cursor with [Shift+R[⌘]] along the time axis, release the mouse button at the end of the section of the mark to be deleted.

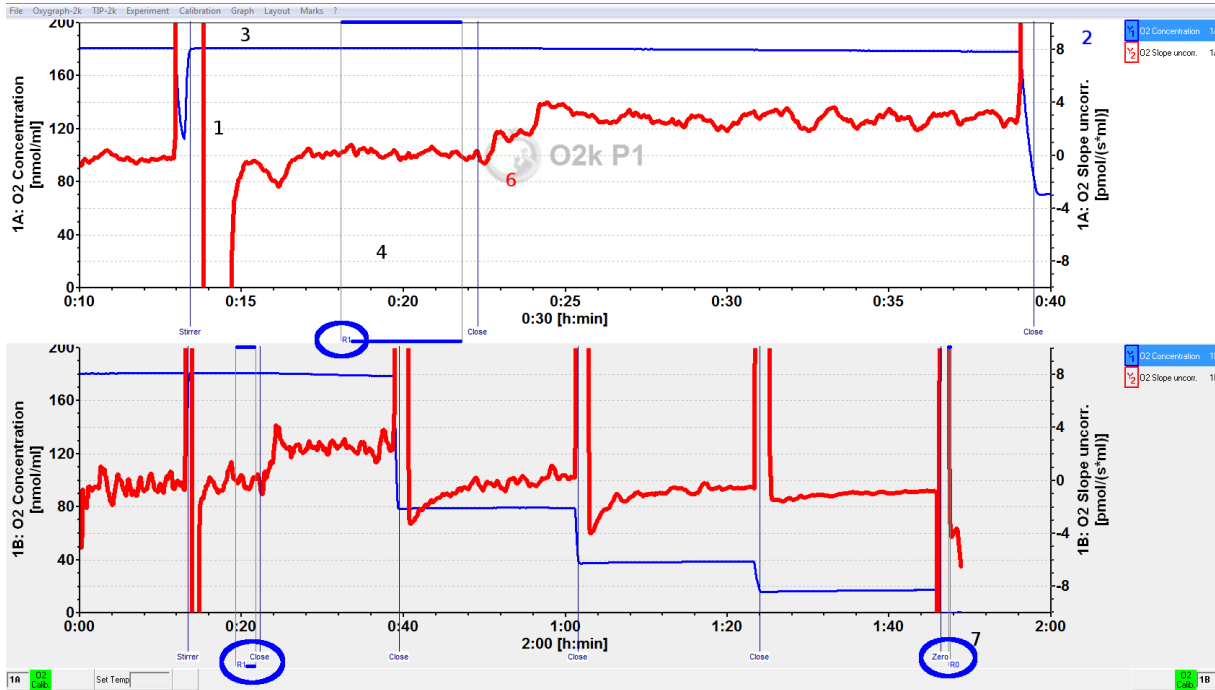

5. Rename the mark: L[Ⓜ] Left mouse click on the bar of the mark. Rename the mark for air calibration as "R1", and the mark for zero calibration as "R0".
6. Observe events (set by [F4]) which indicate particular titrations or any events which are of interest. In the figure, the Event "Close" indicates that the chamber was closed, thus terminating the air calibration phase. Events provide important guidelines for setting and editing calibration marks.
7. Multiple marks can be set.

2.3. DatLab calibration window

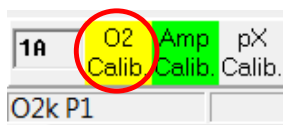

Open the DatLab calibration window: L[Ⓜ] Double click on **O2 Calib.** for O2k-Chamber A or B in the O2k status line.

O2 Calibration

Channel type: 1A: O2 Active POS # 6001

Signal | Slope | Details

Calibration source	Active file	Select Mark	Calib. POS #	POS signal: Recorded [V]	Slope uncorrected [pmol/(s.ml)]	Temperature [°C]	Barometric pressure pb [kPa]
Air calibration: c1	180.97	R1	6001	9.7958	0.17	37.0002	95.20
Zero calibration: c0	0.000	R0	6001	0.0278	0.08	37.0007	95.10

Gain, G [V/μA] 4

O2 solubility factor of medium, FM 0.920 Medium MiR05

Reset to system default

Or open Calibration and select an Oxygen channel. Or select an O₂ Concentration plot and press [F5].

Channel type 1A: O2 Indicates that calibration is performed for the O2 signal in chamber A of Power-O2k P1.

Active POS # The OroboPOS number is displayed as entered in the Oxygraph-2k / O2k Control window ([F7] Tab: Oxygen,

O2) for each chamber. Do not switch sensors between O2k-Chambers, except for troubleshooting.

2.3.1. O2 Calibration \ Tab: Signal

Calibration source The file in which the last calibration has been performed.

Calib. POS # The OroboPOS number for the sensor used in the calibration.

Select Mark ▼ **R1** for "Air calibration": L[⌘] click on the pull down button and select the appropriate mark (R1).

▼ **R0** for "Zero calibration": L[⌘] click on the pull down button and select the appropriate mark (R0). Many times the zero calibration value is used from a previous experiment. The displayed temperature and pressure are without influence on the calibration calculations for zero oxygen.

The average voltage (Raw signal [V]) recorded over the marked section is shown in the corresponding field on the right. The corresponding signal stability is displayed as the uncorrected negative slope of the signal during calibration, Slope uncorr. [$\text{pmol}\cdot\text{s}^{-1}\cdot\text{ml}^{-1}$]. Temperature and barometric pressure are displayed as measured over the marked section.

Calibration values R_1 and R_0 can be edited numerically, without exerting an influence on c_1 . If temperature or barometric pressure are edited, then c_1 is recalculated for the changed conditions.

O₂ solubility factor of medium, F_M Enter the oxygen solubility factor of the medium, F_M , relative to pure water.

More details » [MiPNet06.03 POS-Calibration-SOP](#)


Medium Enter the name of the experimental incubation medium.

c1 The O₂ concentration at air saturation [$\mu\text{mol O}_2\cdot\text{dm}^{-3} = \mu\text{M} = \text{nmol}\cdot\text{cm}^{-3}$], $c_{\text{O}_2}^*$, calculated by DatLab as a function of temperature, barometric pressure, and oxygen solubility factor of the medium.

c0 The known O₂ concentration at the second calibration point. The most practical c_0 is zero oxygen concentration.

Copy from file Calibration parameters can be copied from a file in which a calibration has been performed.

Reset to system default Use only if previously entered and calculated parameters do not make sense.

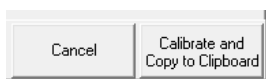

Calibrate and Copy to clipboard After a L[⌘] click on this button, all changes in the entire calibration window

(Tabs Signal and Slope) are applied. The entire plot of oxygen concentration is re-calibrated [$\mu\text{M} = \text{nmol/ml}$], and the corresponding negative slope or volume-specific oxygen flux [$\text{pmol}\cdot\text{s}^{-1}\cdot\text{ml}^{-1}$] is now based on this new calibration. Calibration parameters are automatically copied to clipboard for entry into the "O2k-Calibration-List".

Cancel

Press cancel to exit the calibration window without saving any changes.

2.3.2. O2 Calibration \ Tab: Slope

Slope smoothing ▼ Select the number of data points ($N = 40$ to 5 in intervals of 5) used to calculate the slope through a polynomial fit, as a basis for the plot of flux. A high value of N yields a highly smoothed curve, whereas a low value of N improves time resolution. Apply with **Recalculate slope** or **Calibrate and Copy to Clipboard**. It may take time to recalculate the slope.

Slope amplification Determines the multiplication factor that is applied to the time derivative of the O2 signal.

Background correction Edit the parameters for the background correction of oxygen flux, based on a linear relation between background oxygen flux and oxygen concentration.

a° Intercept [$\text{pmol}\cdot\text{s}^{-1}\cdot\text{ml}^{-1}$]. The default is -2.00 .

b° Slope. The default is 0.025 .

In many applications, background correction based on the default parameters provides sufficient accuracy of respiratory oxygen flux.

O2k-SOP

Background correction is applied automatically to the plots for oxygen flux, except when **Slope uncorrected** is selected. Importantly, for air calibration when the chamber is open for equilibration with a gas phase (air) or for zero calibration, the plot "Uncorrected slope" is selected to evaluate signal stability by observing the zero slope of the oxygen signal.

Further details: [MiPNet14.06](#) and [MiPNet19.18E](#).

Calculate If the active file contains an O2k-background test this function calculates the background parameters, see [MiPNet19.18E](#).

Copy from file Background parameters can be copied from another DatLab file.

2.3.3. O2 Calibration \ Tab: Details

All oxygen calibration parameters are displayed as calculated by DatLab and copied to clipboard upon calibration.

2.4. DatLab calibration: real-time vs. disconnected

DatLab uses calibration values applied real-time (connected to the O2k, recording data) as default values for future experiments. When calibration values are edited in the disconnected mode, they apply only to the current file and will not be used as a new default in experiments. This allows to re-calibrate old files without overwriting the current default values for calibration. Ideally, calibration values that should be used as new defaults are applied real-time when the experiment is still running. However, if the DatLab-calibration is performed after disconnecting, these calibration parameters can be read into other DatLab files using the **Copy from file** function and **Calibrate and Copy to Clipboard**.

Before disconnecting the Oxygraph-2k from DatLab, calibration information is automatically saved and available upon connecting the Oxygraph-2k, even if you exit DatLab and start the program again. The current calibration parameters are displayed when opening the O2 calibration window [F5].



go Bioblast » [MiPNet06.03 POS-Calibration-SOP](#)

3. MultiSensor calibration

Select a MultiSensor plot and press [F5] to open the calibration window, or select the channel from the Calibration menu, or press the Amp or pX buttons in the O2k status line.

pX Calibration		
Channel type:	1A: pX	Active Electrode: 867 Ref: 569
Channel label:	pH	Unit: pH
Signal Slope Details		

Select a Channel and edit **Channel label** and **Unit** If the unit is changed DatLab will ask whether the calibration values should be converted accordingly (Amp calibration only).

Calibrate and Copy to clipboard Apply changes in all tabs (Signal and Slope).



Cancel Press cancel to exit the calibration window without saving any changes.

3.1. Tab: Signal

Calculation of the calibrated signal from the raw signal.

3.1.1. Amp channel: multiple point calibration

The screenshot shows the 'Calibrate' window with the following data:

Select marks	Name	Time	Signal [V]	Slope	Conc. [μM]
<input checked="" type="checkbox"/> 0	0	00:08:58	0.22064	-0.03214	0.00000
<input checked="" type="checkbox"/> 0.1	0.1	00:11:16	0.45685	0.00079	0.10000
<input checked="" type="checkbox"/> 0.2	0.2	00:14:31	0.69102	0.01751	0.20000
<input type="checkbox"/> 0.3					
<input type="checkbox"/> 0					
<input type="checkbox"/> 0.1					
<input type="checkbox"/> 0.3					

Below the table, the 'Gain [V/ μA]' is set to 1000 and 'Polarization voltage [mV]' is set to 100. The 'Sensitivity [V/ μM]' is 2.3519 and 'Intercept [V]' is 0.2210. The 'R²' value is 0.99999. Buttons for 'Copy from file', 'Reset to system default', and 'Show graph' are visible.

In the left panel the current calibration is shown. The right panel is used for a new calibration. Select the marks to use for the calibration (1). The column "Conc." will be filled automatically with the "values" associated with each mark, see [MiPNet19.18](#). Alternatively, enter the known concentrations manually (2a) or

read concentrations from a previous calibration (same mark names) using **Copy from file** (2b). The values associated with the mark will be updated if the concentration was changed in the calibration window. Leave the check boxes beside the Slope column unchecked unless calibrating an Amplex Red experiment:



go Bioblast » http://www.bioblast.at/index.php/Amplex_red

DatLab performs a new linear regression after any field has been edited and displays the sensitivity and intercept (3a). Calibration parameters may be entered manually (3a), or be read from a previous calibration with **Copy from file** (3b), or be **Reset to system default** (3c).

Show Graph Check the quality of the calibration. The graph can be copied to clipboard.

3.1.2. pX channel

Two-point calibration (MiPNet08.16) Linear calibration of pH as a function of recorded voltage is performed by a two-point calibration, using two pH calibration buffers, pX_0 and pX_1 (for example, pH_0 is 7.0 and pH_1 is 4.0; consider the actual pH of calibration buffers at experimental temperature).

Enter pH values or logarithms of two known concentrations in the column labeled "Define two-point

calibration values". Select the corresponding two marks in the "Select Mark" column.

The calibration factor, F , is

$$\text{Eq. 1.} \quad F = \frac{pX_1 - pX_0}{R_1 - R_0}$$

The offset, d , is

$$\text{Eq. 2.} \quad d = \frac{pX_0 \cdot R_1 - pX_1 \cdot R_0}{R_1 - R_0}$$

Calibration of the recorded signal at any time t , R_t , then uses the relation

$$\text{Eq. 3.} \quad pX(t) = (R_t - d) \cdot F$$

Multiple-point calibration

1. Perform a regression of raw voltage [V] against concentration [selected unit] in a spreadsheet program.
2. Note slope and intercept.
3. Open the calibration window.
4. Enter the following data matrix (instead of "Amp conc." "Define two point calibration" shown in the pX calibration window).
5. Press **Calibrate and Copy to clipboard**.



go Bioblast

Calibration of TPP⁺ electrodes » [MiPNet15.03](#).

3.2. Tab: Slope


Controls the calculation of the time derivative (slope) from the signal. In addition to functions described in [Section 2.3](#). [F5] a checkbox allows to calculate slopes from the the raw signal rather than from the calibrated signal.







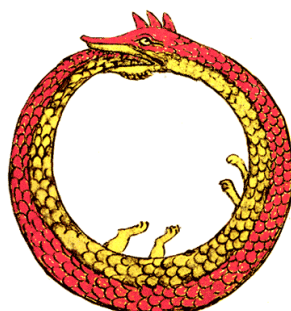

Updates and full version: go Bioblast

» www.bioblast.at/index.php/MiPNet19.18D_O2k-Calibration

Next step – O2k-Manual E » [MiPNet19.18E O2 Flux Analysis](#)



Supplement A. References

- Forstner H, Gnaiger E (1983) Calculation of equilibrium oxygen concentration. In: *Polarographic Oxygen Sensors. Aquatic and Physiological Applications*. Gnaiger E, Forstner H (eds), Springer, Berlin, Heidelberg, New York: 321-3. » 
- Gnaiger E (1983) The twin-flow microrespirometer and simultaneous calorimetry. In: *Polarographic Oxygen Sensors. Aquatic and Physiological Applications*. Gnaiger E, Forstner H (eds), Springer, Berlin, Heidelberg, New York: 134-66. » 
- Gnaiger E (2001) Bioenergetics at low oxygen: dependence of respiration and phosphorylation on oxygen and adenosine diphosphate supply. *Respir Physiol* 128: 277-97. » 
- Gnaiger E (2008) Polarographic oxygen sensors, the oxygraph and high-resolution respirometry to assess mitochondrial function. In: *Mitochondrial Dysfunction in Drug-Induced Toxicity* (Dykens JA, Will Y, eds) John Wiley: 327-52. » 
- Gnaiger E, Steinlechner-Maran R, Méndez G, Eberl T, Margreiter R (1995) Control of mitochondrial and cellular respiration by oxygen. *J Bioenerg Biomembr* 27: 583-96. » 
- Krumschnabel G, Eigentler A, Fasching M, Gnaiger E (2014) Use of safranin for the assessment of mitochondrial membrane potential by high-resolution respirometry and fluorometry. *Methods Enzymol* 542: 163-81.
- Steininger C, Allerberger F, Gnaiger E (2002) Clinical significance of inhibition kinetics in *Streptococcus pyogenes* in response to penicillin. *J Antimicrob Chemother* 50: 517-23. » 



O2k-Manual

- » [MiPNet19.18A](#) Oxygraph-2k: start high-resolution respirometry.
- » [MiPNet19.18E](#) Oxygen flux analysis: DatLab real-time.
- » [MiPNet15.03](#) O2k-MultiSensor system with ion selective electrodes (ISE).
- » [MiPNet15.05](#) O2k-Manual: amperometric sensors (NO).



O2k-Protocols: O2k-SOPs

- » [MiPNet06.03](#) POS calibration SOP.
- » [MiPNet14.06](#) Instrumental background and accuracy of oxygen flux.

Supplement B. O2 Calibration with titanium stoppers



We recommend to replace titanium stoppers with the new PEEK stoppers and corresponding O2k-Chamber holders.

Stopper-spacers used for air calibration with the PVDF stopper (Chamber A) or titanium stopper (Chamber B).

O2k-Catalogue:

- » http://wiki.orooboros.at/index.php/O2k-Chamber_Holder
- » http://wiki.orooboros.at/index.php/Stopper%5Cblack_PEEK%5Cconical_Shaft%5Ccentral_Port