

# Data analysis of mitochondrial membrane potential estimation using various fluorescence dyes

Komlodi T<sup>1</sup>, Antunes D<sup>1</sup>, Iglesias-Gonzalez J<sup>1</sup>, Sobotka O<sup>2,3</sup>, Gnaiger E<sup>1</sup>

## <sup>1</sup> Oroboros Instruments

Schoepfstrasse 18, A-6020 Innsbruck, Austria

Email: [instruments@oroboros.at](mailto:instruments@oroboros.at)

[www.oroboros.at](http://www.oroboros.at)

<sup>2</sup> 3rd Department of Internal Medicine, Gerontology and Metabolism  
University hospital and Faculty of Medicine in Hradec Kralove  
Charles University, CZ

<sup>3</sup> Department of Physiology, Faculty of Medicine in Hradec Kralove  
Charles University, CZ

## Contents

1. General information .....	1
2. Starting data analysis .....	2
3. Calibration .....	3
4. Chemical background .....	3
5. Biological Experiment .....	3
5.1. Oxygen flux analysis .....	3
5.2. Membrane potential analysis .....	4
6. References .....	5

## 1. General information

**Substrate-uncoupler-inhibitor-titration (SUIT) protocols** are designed to study respiratory control in a sequence of coupling and pathway control states induced by multiple titrations within a single experimental assay. DatLab 7.4 has been specifically designed to guide the user through SUIT protocols ([DL-Protocols](#) in DatLab). Excel templates (SUIT-###\_Fluo\_mt\_D###\_general.xlsx) are provided for data analysis of O<sub>2</sub> flux and relative mitochondrial membrane potential (mtMP) using different fluorescence dyes (*e.g.*, safranin, TMRM, Rhodamine123) for isolated mitochondria, tissue homogenate and permeabilized cells. Each DL-Protocol is defined with a unique D-number (D###), for a detailed list see:

- [https://www.bioblast.at/index.php/SUIT\\_protocol\\_library#List of S  
UIT protocols with D-numbers](https://www.bioblast.at/index.php/SUIT_protocol_library#List_of_SUIT_protocols_with_D-numbers)


Use the SUITbrowser to find the best SUIT protocol for your research questions:

- <https://suitbrowser.oroboros.at/>

## 2. Starting data analysis

Upon completion of real time respirometry measurement in DatLab 7.4 or anytime while doing the DatLab analysis, open our Excel template to analyze your data in a time efficient way.

1. In DatLab 7.4, select the menu [Protocols] and click on [SUIT:Browse DL-Protocols and templates].

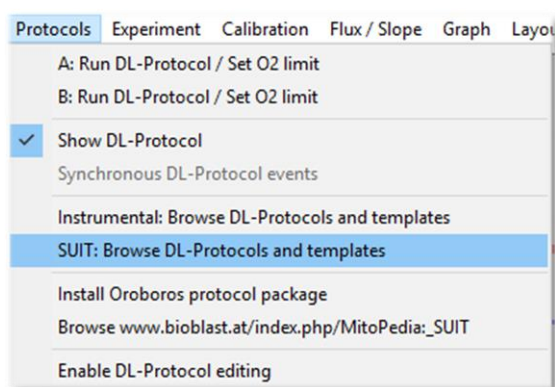

2. Select your SUIT protocol and open the SUIT-###\_Fluo folder. Inside this folder, you will find another folder for the specific DL-Protocol (named SUIT-###\_Fluo\_mt\_D###). In each folder, four Excel files can be found:
  - a. A blank template (named SUIT-###\_Fluo\_mt\_D###\_general.xlsx) to calculate the relative mtMP values.
  - b. A demo version of the general template (named SUIT-###\_Fluo\_mt\_D###\_general\_demo.xlsx), which provides an example of the file already with data showing the relative mtMP values.
  - c. A blank template (named SUIT-###\_Fluo\_mt\_D###\_safranin.xlsx) to convert the fluorescence values measured by safranin into absolute mtMP values expressed as mV.
  - d. A demo version of the specific safranin template (named SUIT-###\_Fluo\_mt\_D###\_general\_demo.xlsx) which provides an example of the file already with data showing the mtMP values expressed in mV.

3. Create a copy of SUIT-###\_Fluo\_mt\_D###\_general.xlsx analysis template for your data analysis and rename it. You can rename the template by opening it and choosing the option 'Save as' in the archive top menu.

### 3. Calibration

- 3.1. Open the DatLab file containing the data from the calibration.
- 3.2. Open the menu [Calibration] and select 'Amperometric, Amp' to calibrate and convert the amperometric signal into the concentration of the fluorescence dye.
- 3.3. Select the first four marks (e.g., Saf0, Saf0.5, Saf1, Saf1.5, Saf2) for calibration. Check the  $r^2$  of the linear regression and press the 'Show graph' button to check the linearity of regression. The sensitivity [ $V/\mu\text{M}$ ] can be also found in the same window.
- 3.4. Press 'Calibrate' and the [Y1: Amp raw] will turn into calibrated [Y1: Amp]. Adjust the scaling [F6] of [Y1: Amp].

### 4. Chemical background

Please follow instructions in section 4 of [MiPNet24.08 Safranin Analysis Template](#).

### 5. Biological Experiment


In the Excel template you can select the setting by ticking the boxes 'Titration volume correction' and 'Known sample concentration'. More information can be found here: [MiPNet24.06 Oxygen flux analysis with DatLab7.4](#)

#### 5.1. Oxygen flux analysis

The calculations of the  $\text{O}_2$  fluxes are provided under the following link complying with Oroboros transparency policy:

<https://wiki.orooboros.at/index.php/Flux / Slope#O2>

- 4.1.1. In DatLab 7.4, after setting the marks separately to the  $\text{O}_2$  flux, go to [Marks], and select 'Slope uncorrected + all info' in 'DL-Protocol marks'. In the new window select:
  - a. Your chamber of interest.
  - b. Plot for Marks: 'O2 slope neg. [ $\text{pmol/s}\cdot\text{mL}$ ]'.
  - c. Channel: 'Oxygen,  $\text{O}_2$ '. Leave only this channel selected.
  - d. Select: 'Median'.
  - e. Sort by: 'Time' (default).
  - f. Then, click on [Copy to clipboard] to copy the selected values.

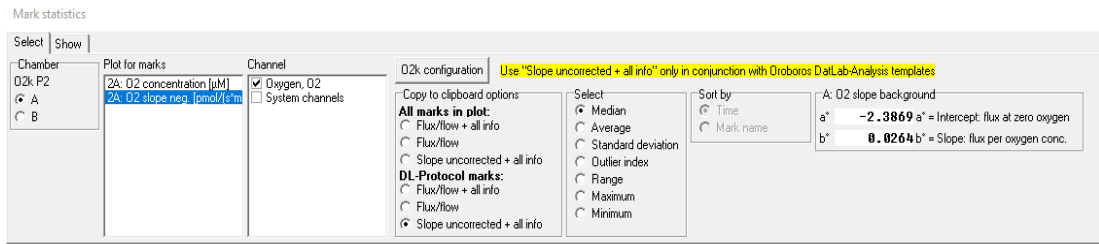

- 4.1.2. In the Excel template (SUIT-###\_Fluo\_mt\_D###\_general.xlsx): Click on the yellow cell A7 and paste these selected values (only O<sub>2</sub>) from DatLab [Ctrl+V].
- 4.1.3. The calculated values for the specific O<sub>2</sub> flux, specific O<sub>2</sub> flux (bc), *FCR* and *FCR* (bc), on each step of the protocol can be found from column K in the rows 24, 25, 27 and 28, respectively.
- 4.1.4. Paste the DatLab graphs showing the traces for the chamber:
- In DatLab: Select the graph (left mouse click into the respirometry graph of interest) → select 'Graph\Copy to Clipboard\WMF'.
  - In the Excel template: Click on the yellow cell A27: 'Paste DatLab graph here, reduce to width 22 cm (8 inches)' → press [Ctrl+V] to paste.
  - Select the graph (right click on the graph) → select 'Size and properties' and set the width of the graph to 22 cm (8 inches).

## 5.2. Membrane potential analysis

- 5.2.1. In DatLab 7.4, copy the calibration values of the fluorescence signal from the previous calibration file: Open the menu [Calibration] and select 'Amperometric, Amp', press 'Copy from file' to open the calibration file and copy the sensitivity value. Then press the 'Calibrate' button in the 'Amp calibration' window.
- 5.2.2. Select [Y1: Amp] as the active plot for setting marks and place marks according to your protocol as done for the O<sub>2</sub> flux.
- 5.2.3. Open [Marks] window, select 'DL-Protocol marks' and 'Slope uncorrected + all info'; in the new window select channel 'Amperometric, Amp' and use DL-Protocol marks: 'Slope uncorrected + all info' to display data for the marked regions.
- 5.2.4. Export data with 'Copy to Clipboard'.
- 5.2.5. Paste the calculated fluorescence values in the Excel template into the yellow cell A42.
- 5.2.6. The calculated relative mtMP values can be found in row 62 starting with column L.
- 5.2.7. Copy the Amp graph as it is explained in the 4.1.4 and paste on the yellow cell A63.

## 6. References

1. Cardoso LHD, Antunes D, Iglesias-Gonzalez J, Komlodi T, Doerrier C, Garcia-Souza LF, Gnaiger E, Sobotka O (2019) Oxygen flux analysis with DatLab 7.4. Mitochondr Physiol Network 24.06(01):1-5. - [»Bioblast link«](#)
2. Krumschnabel G, Fasching M, Gnaiger E (2019) O2k-FluoRespirometry: HRR and simultaneous determination of mt-membrane potential with safranin or TMRM. Mitochondr Physiol Network 20.13(03):1-5. - [»Bioblast link«](#)



» [MitoPedia:\\_Mitochondrial\\_membrane\\_potential](#)

» [Flux / Slope](#)

» [Safranin](#)

» [TMRM](#)

» [Rhodamine123](#)

## Acknowledgements



The project NextGen-O2k has received funding from the European Union's Horizon 2020 research and innovation programme under grant agreement No 859770.

